

StarLab Release Notes

Version 4.00 (.08) Released 12-May-2026

New features:

- Support for automatic conversion of Ariel logfiles uploaded by StarLab or StarViewer, and native logfiles from StarViewer. Reads and displays the following log file types:
 - StarLab native log
 - StarLab-uploaded Ariel log
 - StarViewer/Android native log
 - StarViewer/Android-uploaded Ariel log
 - StarViewer/iOS-uploaded Ariel log

Updates:

- Updated "Ophir User Commands"
- Updated build files to use Visual Studio 2022
- MSVC v14 Redistributable that comes with VS 2022
- Minor internal updates

Fixes:

- Log upload progress bar fixed
- Minor internal fixes

Known Issues:

- Time-out when searching for non-existing devices on Bluetooth or Ethernet can take up to a minute!
- COM-Object bug mainly affecting Python applications

Version 3.93 (.02) Released 23-December-2024

New features:

- The COM object now supports wireless connections to Ariel

Updates:

- COM objects version 10.8
- Updated "Ophir User Commands"

Fixes:

- Allow DOSAGE mode for StarLite, with PD300-RM

Known Issues:

- Time-out when searching for non-existing device on Bluetooth or Ethernet can take up to a minute!

Version 3.92 (.03) Released 11-July-2024

New features:

- Added PD300RM support for Centauri on StarLab
- Enhancement: Add auto scale option for line graphs
- Support connecting to Centauri via Ethernet
- Extended "Time Frame" of graphs to be up to 99:59 hours

Updates:

- Update "Ophir User Commands"
- COM objects version 10.7

Fixes:

- StarLab previous version didn't run on Win 7
- When attempting to open a second instance of StarLab there was an assertion failure

Known Issues:

- Time-out when searching for non-existing device on Bluetooth or Ethernet can take up to a minute!